SONOPULS

Ultraschall-Homogenisatoren

Ultrasonic Homogenizers

Emulgieren

Homogenisieren

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Suspendieren

Sonochemie

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Emulsifying

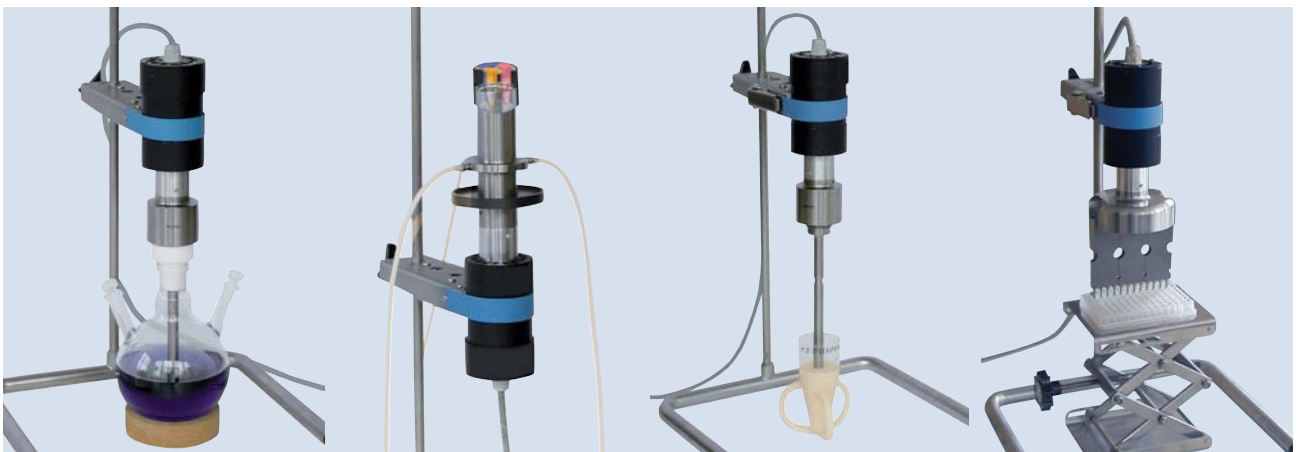
Homogenizing

Cell Disruption

Suspending

Sonochemistry

Accelerating Reactions

BANDELIN

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55 Years of Experience in Ultrasound Technology

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How to select the proper unit

Power output in watt is not the sole criterion for selecting an ultrasonic homogenizer. This value only indicates the power of the HF-generator but not the energy delivered to the sample. The amplitude at the radiating surface of the probe is the determining factor for the evaluation of the irradiation result while taking into consideration the volume of the sample.

SONOPULS ultrasonic homogenizers produce with the same electric power higher amplitudes than other models in the market thanks to an ideal integration of all components.

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AMPLICHRON®-system guarantees a constant amplitude independently from changing conditions within the sample. It ensures reproducible results for process validation. Settings within a range of 10 to 100 %. Verification of actual value at the display. Permanent control of ultrasound irradiation as well as indication of wear of the probe.

Pulsation

limits increase of temperature when processing sensitive samples. The adjustable pulsation allows cooling during rest intervals.

Built-in timer

Process duration storable. Indication of elapsed time in continuous operation or of remaining time in countdown mode.

Processing frequency - 20 kHz uniform sonic waves

Constant sound radiation- extremely effective.

Switching ON / OFF - easy to handle

either at the generator or directly at the ultrasonic converter via button or remote control.

Accessories

A wide range of probes and special accessories for a vast variety of applications

Foil keypad

easy to clean and user-friendly

Fail-safe during continuous operation and idling

RFI-proofed and CE-marked, also as medical device compliant to the directive for in-vitro diagnostics 98/79/EG

Mode of operation

HF generator:

Transforming of low-frequency voltage of 50 Hz into high-frequency voltage of 20 kHz.

Ultrasonic converter:

Transforming of electrical energy delivered from the generator into mechanical vibrations of 20 kHz.

Standard and booster horns:

Horns increase through their specially designed shape the amplitude. The external thread is made for close connection of vessels.

Probes:

Transmitting of ultrasonic energy into the sample. Microtips, tapered and flat tips, dia. 2, 3, 6, 13, 19 and 25 mm, for use in different volumes.

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Small unit for lab routine

Ready-to-operate basic equipment for volumes from 2 ml to 50 ml consisting of generator GM 2070, ultrasonic converter UW 2070, stepped standard horn SH 70 G and microtip MS 73, 3 mm diameter. HF-output max. 70 W_{eff}

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Standard unit for lab routine

Ready-to-operate basic equipment for volumes from 20 ml to 900 ml consisting of generator GM 2200, ultrasonic converter UW 2200, booster horn SH 213 G and titanium flat tip TT 13 of 13 mm diameter. HF-output max. 200 W_{eff}

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High tech for research

- for small volumes -
Ready-to-operate basic equipment for volumes from 2 ml to 50 ml consisting of generator GM 3100, ultrasonic converter UW 3100, stepped standard horn SH 70 G and microtip MS 73, 3 mm diameter. HF-output max. 100 W_{eff}

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High tech for research

- for larger volumes -
Ready-to-operate basic equipment for volumes from 20 ml to 900 ml consisting of generator GM 3200, ultrasonic converter UW 3200, booster horn SH 213 G and titanium flat tip TT 13 of 13 mm diameter. HF-output max. 200 W_{eff}

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High tech for research and pilot plant stations

Ready-to-operate basic equipment for volumes from 100 ml to 2500 ml consisting of generator GM 3400, ultrasonic converter UW 3400, booster horn SH 3425 and extended probe, 25 mm diameter. HF-output max. 400 W_{eff}

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Waste water samples

Aim: Homogenizing for determination of harmful substances, e.g. mineral oil, grease AOX in industrial and butcher's waste water

Quantity: 250 ml

Approx. time: 5 - 10 min

Unit: HD 2200/3200 with TT 13, for taller vessels with VS 70 T

Aluminium oxide suspensions

Aim: Dispersing

Quantity: 100 ml

Approx. time: ca. 4 min

Unit: HD 3200 with KE 76

Soil samples

Aim: Extraction for determination of pH value, Mg, K, P, N – contents for recommendation of fertilizer / determination of radio nucleides to control radioactivity in the environment (milk researchment)

Quantity: 50-100 ml / 100-150 ml

Approx. time: a few seconds

Unit: HD 2200/3200 with KE 76 / VS 70 T

Bladder tissue

Aim: Disruption

Quantity: 1,5 ml

Approx. time: ca. 1,5 min

Unit: HD 2200 with MS 72, cooling necessary

Candida albicans

Aim: Disruption

Quantity: 10 ml

Approx. time: 10 min

Unit: HD 2070/3100 with MS 73

Large intestine tissue

Aim: Disruption

Quantity: 1,5 ml

Approx. time: ca. 3 min

Unit: HD 2200 with MS 72, cooling necessary

Dispersing of solid particles

Aim: grain size analysis

Quantity: 50 - 100 ml

Approx. time: 2 - 5 min

Unit: HD 2200/3200 with KE 76

Small intestine tissue

Aim: Disruption

Quantity: 1,5 ml

Approx. time: ca. 2,5 min

Unit: HD 2200 with MS 72, cooling necessary

Escherichia coli

Aim: Disruption for proteine lay off

Quantity: 10 ml

Time: 5 - 10 min

Unit: HD 2070/3100 with MS 73 or HD 2200 with MS 73

Eucaryotic cells

Aim: Disruption for proteine lay off

Quantity: 1,5 ml

Approx. time: 1 min

Unit: HD 2200/3200 with BR 30 and EH 3, cooling necessary

Meat and sausage samples

Aim: Homogenizing for determination of nitrates

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Insect cells

Aim: Disruption for proteins lay off
Quantity: 20 - 50 ml
Approx. time: app. 25 sec, pulsed
Unit: HD 2070/3100 with MS 73 and RZ 2

Liver tissue

Aim: Homogenizing for molecular genetic tests
Quantity: 1,5 ml
Approx. time: 1½ min
Unit: HD 2200/3200 with MS 72, cooling necessary

Liposomes

Aim: Producing of small unilamellar phospholipid vesicles
Quantity: 20 ml
Approx. time: 10 - 15 min
Unit: HD 2070/3100 with TT 13, cooling necessary

Lymphocytes

Aim: Disruption
Quantity: 50 µl - 2 ml
Approx. time: 1 - 5 min
Unit: HD 2070/3100 with BR 30 and EH 3

Nano emulsions

Aim: Drop sizes within nm range
Quantity: 2 ml
Approx. time: ca. 5 min
Unit: HD 3100 with MS 72, cooling necessary

Nano particles

Aim: Dispersing
Quantity: 1,5 ml
Approx. time: 4 min
Unit: HD 3200 with KE 76

Kidney tissue

Aim: Homogenizing
Quantity: 1,5 ml
Approx. time: ca. 40 sec
Unit: HD 2200/3200 with MS 72, cooling necessary

O/W emulsions

Aim: finest emulsions
Quantity: 10 ml
Approx. time: ca. 1 min
Unit: HD 3200 with KE 76,
vessel: rosett cell

Carbon black dispersions

Aim: Homogenizing
Quantity: 50 ml
Approx. time: ca. 5 min
Unit: HD 2200 with DH 13 G, vessel: KG 3

Saccharomyces cerevisiae

Aim: Disruption
Quantity: 20 ml
Approx. time: 15 min
Unit: HD 2200/3200 with KE 76, addition of glass beads to accelerate process, cooling necessary

Staphylococcus aureus

Aim: Disruption
Quantity: 10 ml
Approx. time: 10 min
Unit: HD 2070/3100 with MS 73

Streptococcus

Aim: Disruption
Quantity: 10 ml
Approx. time: 8 - 10 min
Unit: HD 2200/3200 with MS 73

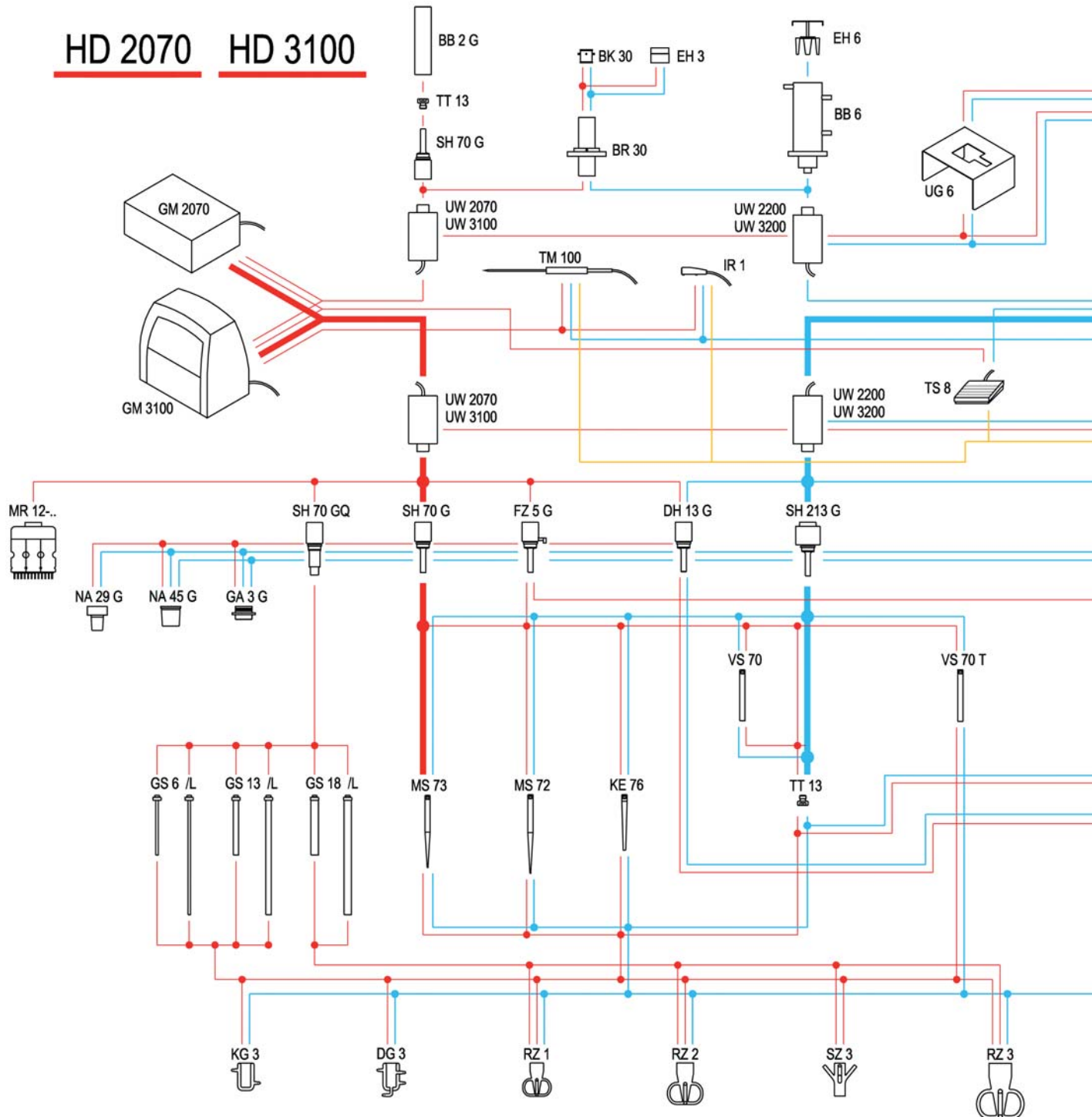
Detection of prions through cyclic amplification of protein misfolding

Quantity: 200 ml
Approx. time: ca. 1,5 min
Unit: HD 2070 with MS 73

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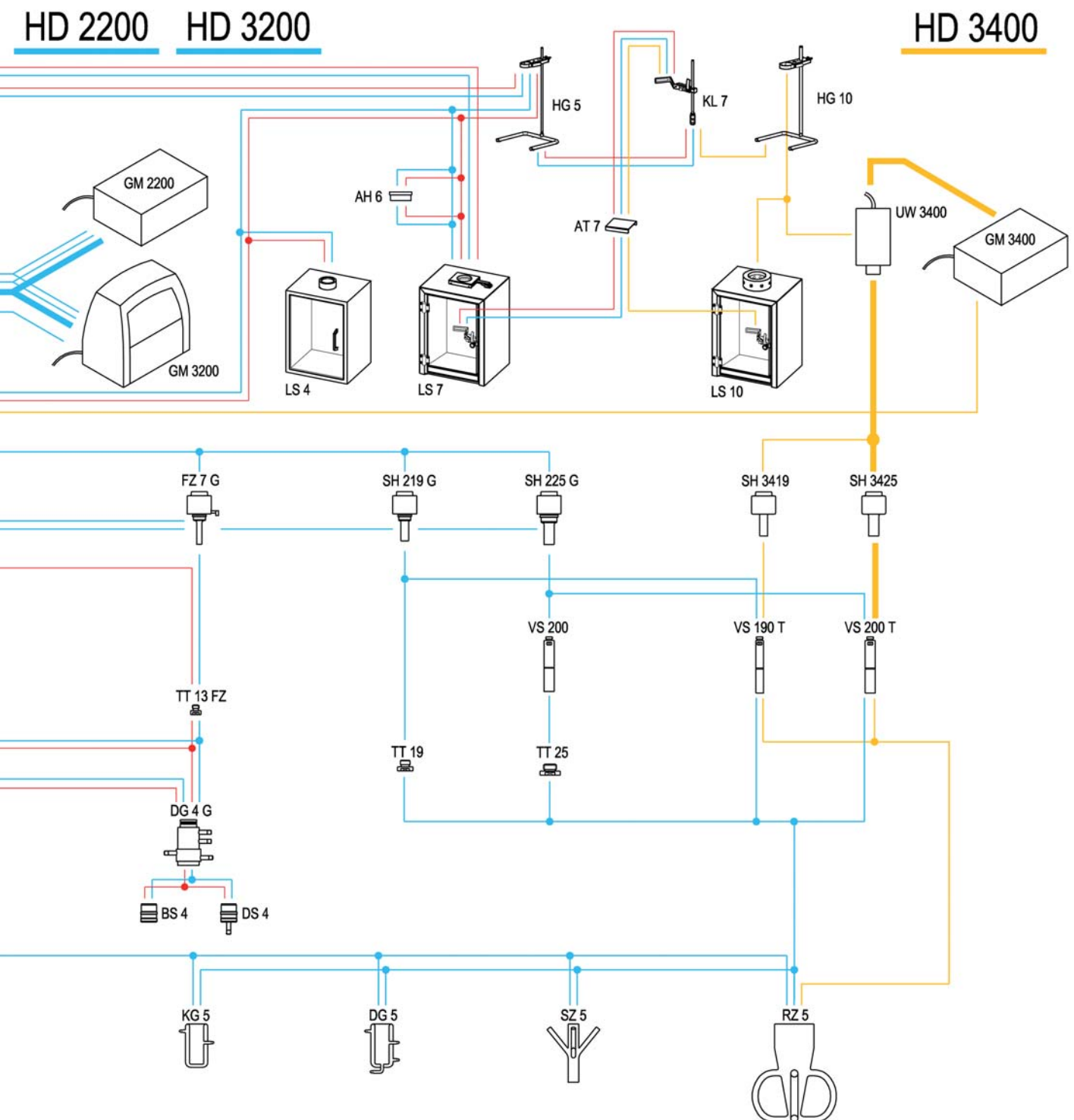
Accessories and Applications

Ultrasonic homogenizers are used in laboratories, hospitals and in industry for scientific experiments and analysis as well as in pilot or small-lot production. Here are some examples showing the vast variety of applications for ultrasonic homogenizers:

- Disruption of cells, bacteria, viruses, tissue
- Emulsifying, e. g. in cosmetics
- Homogenizing of samples
- Preparing of samples for HPLC
- Degassing of liquids
- Preparing of test portions for determination of grain size
- Acceleration of chemical reactions
- Analysis of waste water

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Probes made of titanium alloy (Ti-Al6-V4) transmit mechanical longitudinal waves into the sample. They are thermo-resistant, can be treated in autoclaves and are resistant to corrosive media. Sample volume, diameter of the processing vessel and the required amplitude determine the selection of the unit and the type of probe. The higher the amplitude, the more intense the sonication. For sonication of suspensions only long probes should be used. Safe application – no suspension particles in the thread.

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VS for enlarging the operating depth when using flat tips.
VS 70 between SH 70 G / 213 G and TT 13
VS 200 between SH 225 G and TT 25

Typ / type	VS 70	VS 200
für / for HD	2070/2200/3100/3200	2200/3200/3400
Best.-Nr. / Code No.	500	415

Bezeichnung description		Mikrospitzen microtips		Kegelspitze tapered tip	Sonotrode lang extended probe		
Typ / type		MS 72	MS 73	KE 76	VS 70 T	VS 190 T	VS 200 T
Best.-Nr. / Code No.		492	529	530	494	3638	478
Durchmesser / diameter	mm	2	3	6	13	19	25
Länge ca. / length approx	mm	191	175	135	126	121	139
Stufenhorn für / standard horn for HD 2070/3100		SH 70 G	SH 70 G	SH 70 G	SH 70 G	-	-
Boosterhorn für / booster horn for HD 2200/3200		SH 213 G	SH 213 G	SH 213 G	SH 213 G	SH 219 G	SH 225 G
Boosterhorn für / booster horn for HD 3400		-	-	-	-	SH 3419	SH 3425
Amplitude für / amplitude for HD 2070/3100	µm _{SS} *	253 / 285	212 / 245	165 / 191	80 / 97	-	- / -
Amplitude für / amplitude for HD 2200/3200		282 / 286	302 / 308	249 / 255	153 / 170	73 / 81	46 / 51
Amplitude für / amplitude for HD 3400		-	-	-	-	116	82
Volumen / volume HD 2070 / 3100	ml	1 - 25	2 - 50	5 - 100	10 - 200	-	-
Volumen / volume HD 2200 / 3200	ml	2 - 30	5 - 90	10 - 350	20 - 900	25 - 900	30 - 1000
Volumen / volume HD 3400						500 - 1500	500 - 2500
Gefäß Ø mind. / vessel diameter min	mm	4	6	8	17	23	29

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Standard and booster horns (Ti-Al6-V4) are furnished with a thread for replaceable probes. With exterior thread to connect various vessels.

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Solid standard horn
DH 13 G with diamond coating on the radiating area; lifetime is thirty times longer than usually.

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MR for connection to HD 2070/3100.
*Composed of Multison horn MRH 12
and 12 Multison tips MRS 2, MRS 3 or
MRS 2-2C .*
*For irradiation of samples in microplates and
deep well plates.*
Simultaneous sonication of 12 samples.
Multison tips individually replaceable .

Typ / type	MR 12-2	MR 12-3	MRH 12-2C
Best.-Nr. / Code No.	3626	3633	3643

Typ / type	MRS 2	MRS 3	MRS-2C
Ø / dia.	2 mm	3 mm	2 mm
Material / material	Ti6Al4V	Ti6Al4V	Quarzglas / silica glass
Best.-Nr. / Code No.	3628	3629	3642

Typ / type	MRH 12
Best.-Nr. / Code No.	3627

Titanteller <i>titanium flat tips</i>			Quarzglas-Sonotroden <i>silica glass probes</i>					
TT 13	TT 19	TT 25	GS 6	GS 6 L	GS 13	GS 13 L	GS 18	GS 18 L
497	491	532	024	048	028	050	040	054
13	19	25	6		13		18	
5	5	6	145	290	145	290	145	290
SH 70 G	-	-	SH 70 GQ		SH 70 GQ		SH 70 GQ	
SH 213 G	SH 219 G	SH 225 G	-		-		-	
-	-	-	-		-		-	
78 / 93	- / -	- / -	12,5 / 13		13 / 13		13 / 13	
149 / 165	73 / 81	48 / 53	- / -		- / -		- / -	
-	-	-	- / -		- / -		- / -	
10 - 200	-	-	2 - 100		25 - 200		25 - 500	
20 - 900	25 - 900	30 - 1000	-		-		-	
-	-	-	-		-		-	
17	23	29	10		17		22	

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GS for connection to HD 2070/3100 with
special horn SH 70 GQ.
*For application in food analysis, pharmacy or
environmental analysis. No intrusion of metal
particles and boron compounds - ideal for
trace analysis. High chemical and temperature
shock resistance, no electric conductivity.*

Typ / type	SH 70 GQ
Best.-Nr. / Code No.	013

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FZ 5 G and FZ 7 G to prepare stable mixtures
of non-mixable or hardly mixable liquids (oil-in-
water) by direct intrusion of pre-mixed samples
into the cavitation field. In combination with
flow-through cell DG 4 G the continuous treat-
ment of 2 different media and parallel tem-
pering is possible.
Material: Ti-6Al-4V

Typ / type	FZ 5 G	FZ 7 G	TT 13 FZ
für / for HD	incl. TT 13 FZ	incl. TT 13 FZ	für/for FZ 5 G / FZ 7 G
Best.-Nr. / Code No.	2070/3100 490	2200/3200 452	496

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Processing vessel, stainless steel
DG 4 G for high-volume flow-through processing like emulsifying, dispersing or homogenizing, up to 30 l/h,
KG 4 G, closed vessel with cooling jacket. Processing volume about 65 ml.

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Processing vessels made of glass
Cooling vessel KG for sonication of temperature-sensitive samples.
Flow-through vessel DG with cooling jacket for irradiation of larger volumes.
Rosett cell RZ for homogenous and intense circulation of sample.
Suslick cell SZ with 3 sidearms for introduction of gas or measuring probes.

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Sleeve adapters made of PTFE for tight mounting on standard ground glass vessels.
NA 29 G for NS 29/32 for SH 70/213 G
NA 45 G for NS 45/40 for SH 70/213/219/225 G
GA 3 G threaded adapter made of stainless steel with external thread M 40 x 1 for SH 70/213/219/225 G

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Microtube holder EH 6

For use in BB 6. Up to 6 samples can be treated simultaneously. Pressure plate holds tubes in place.
No floating of cups.

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Cup Horn BB 6

Temperature regulation with circulating liquid coolants

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Microtube holder EH 3 for use with BR 30. Up to 3 samples can be treated simultaneously. 2 exchangeable discs with diameters 8,5 or 11,5 mm.

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Inset basket BK 30

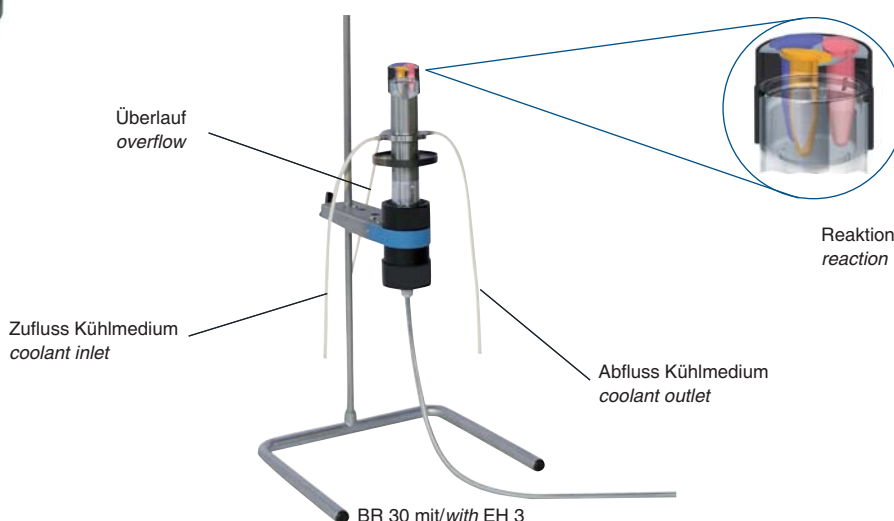
For intensive cleaning of small parts, e. g. cleaning of radioactively contaminated seeds in BR 30.

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Cup booster BR 30

For high-intensive irradiation of smallest and sensitive sample volumes.

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Typ / type HG 5
für / for HD 2070/2200/3100/3200
Best.-Nr. / Code No. 459

Stainless steel stand
with lab clamp, for safe fixing of ultrasonic converter.

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Best.-Nr. / Code No. 3636

Clamping device KL 7 (DE 20 2006 005 654.98) for HG 5 / HG 10 with rod, swivelling clamp and quick clamp for reaction vessels dia. 15 mm to dia. 100 mm

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Best.-Nr. / Code No. 3644

Supporting table AT 7 suitable for KL 7 or in LS 7 with non-slip mat to hold sample vessels safely in place

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LS 4
kunststoffbeschichtete Spanplatten, 10 dB-AU Dämpfung.

Sound proof boxes reduce the noise level considerably. Precut holes at the backside allow connections for gas supply and flow-through processing. Acrylic door permits process monitoring.

LS 4
Plastics coated walls, 10 dB-AU damping.

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LS 7
Made of stainless steel, with damping plates. 20 dB-AU damping.

With rod, swivelling clamp and quick clamp for height adjustment of sample vessels.

Clamping belt for safe fixing of sample vessels with different sizes. Also applicable for sonication of samples in glass vessels with round bottoms or with inlets from below.

Special support UG 6 is available for inverted position of the box with cup horn BB 6 or cup booster BR 30. Ultrasonic converter is fixed safely through a special clamp.

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LS 10
Made of stainless steel, with damping plates. 20 dB-AU damping.

LS 4	LS 7	LS 10	UG 6
2070/2200	2070/2200	3400	2070/2200
3100/3200	3100/3200		3100/3200
10	20	20	
416	3635	3637	3616

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Typ / type AH 6
für / for LS 7
Best.-Nr. / code no. 3619

Distance tube for direct sonication with long probes.

AH 6: MS 72/73, KE 76, VS 70 with TT 13, VS 200 with TT 25 / VS 200 T, VS 70 T, GS ...

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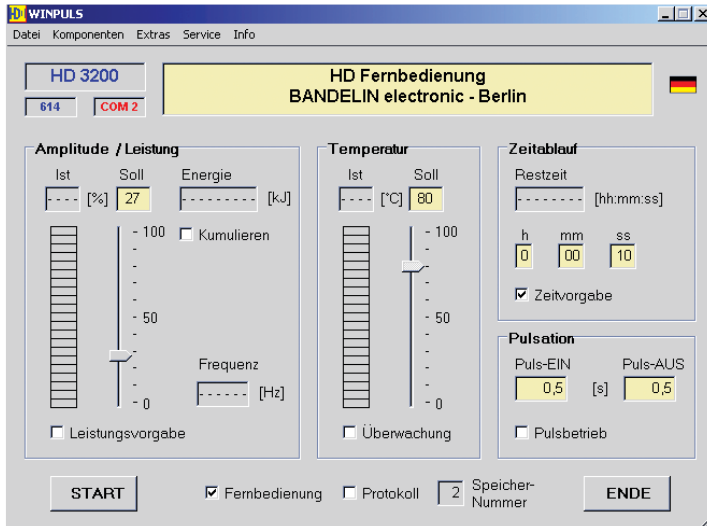
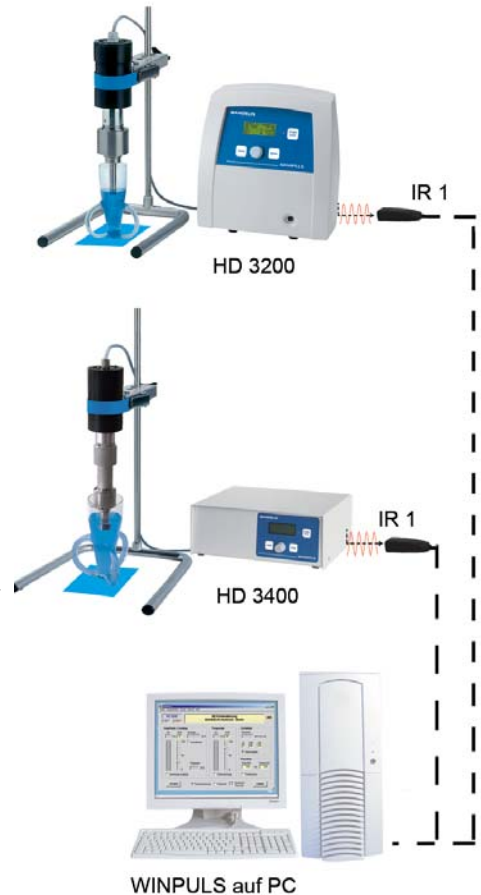
For process control with PC with operating systems WINDOWS® 2000 and WINDOWS® XP. With different additional functions like test logging and comfortable data storage (up to 99 storages). Set composed of WINPULS® software and infrared adapter IR 1 for interface RS 232.

Typ / type WINPULS®-Fernsteuerung mit Infrarot-Adapter IR 1
WINPULS® remote control with infrared adapter IR 1
für / for HD 3100/3200/3400
Best.-Nr. / Code No. 3625

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
Typ / type TS 8
für / for HD 2070/2200
3100/3200/3400
Best.-Nr. / Code No. 531

Foot switch remote control for easy switching ON/OFF of the HF-generator. With 3 m cable.

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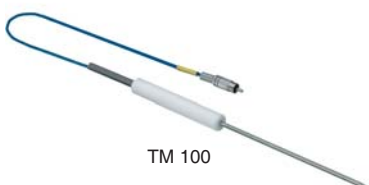Typ / type TM 100
für / for HD 3100/3200/3400
Best.-Nr. / Code No. 3622

Temperature sensor for measuring of sample temperature from 0 up to 120 °C.

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Examples of applications

- *Disinfection of organic substances in rinsing liquids for recycling*
- *Dispersion of nano-scaled polishing suspensions used in wafer industries*
- *Dispersion of nano-porous clay particles in inkjet paper manufacturing*

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*Consisting of:
Cylindrical immersible transducer
RT 4-1040
Reactor housing RG 4-000
HF Generator LG 1001 T, 1000 W*

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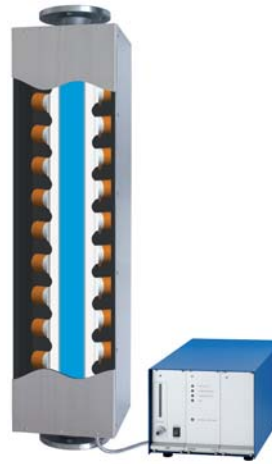
Examples of applications

- *Sewage sludge disintegration*
- *Production of ceramics suspensions*
- *Dispersion of silicic acid used in wafer industries*
- *Production of PTFE suspensions used for coatings*

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- *standard baths from 1 - 200 litres*
- *compact construction*
- *easy to handle*
- *optionally with heating*
- *integrated timer*
- *from 3 ltr. upwards with drain*
- *cost-saving - reduced cleaning time*
- *intensive, gentle, thorough*
- *large range of accessories*

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BANDELIN *electronic*
specialized in manufacturing of ultrasonic units, maintain a quality management system complying with the requirements of
EN ISO 9001:2000
and EN ISO 13485:2003

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GmbH & Co. KG