



Electronic Crimpers and Decappers

Hand-held Rechargeable Crimpers and Decappers

- Precision, rechargeable crimpers and decappers for 8, 11, 13 & 20mm caps
- Easy to use, comfortable grip
- Robust, ergonomic design for left or right handed use
- Hundreds of reproducible crimps from one charge
- Crimp force sensing assures consistent proper sealing
- Adjustable crimp settings for compatibility to most cap types




Reliable, Reproducible Crimping with the Push of a Button

**Tired of crimping hundreds of vials with manual crimpers?
Tired of losing your samples due to poorly crimped seals?**

The motor driven, Electronic Crimper solves these problems and crimps hundreds of vials on a single charge with just the touch of a button.

The electronic Crimpers and Decappers are hand held tools for attaching or removing aluminum seals on 8mm, 11mm, 13mm and 20mm vials. A single press of the button both closes the crimping jaws to attach the seal and opens them to release the vials. You can also remove a seal with a single press of the operation button. The tools are compatible with standard aluminum seals, pressure release seals, tin plated seals and bi-metallic two piece seals. (Select the proper tool according to the finish size of the vials). The long life lithium ion battery provides full power crimping and decapping throughout the charge cycle.

* Flip top seals with plastic buttons on the top are not compatible with the electronic crimpers.

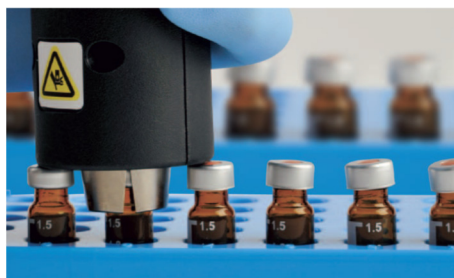

Features

- Secure, reproducible crimps with the push of a button
- Quick and easy seal removal with the push of a button
- Ergonomic push button design eliminates wrist strain
- Long life lithium ion battery with low battery indicator
- Universal 100V-240V charger included
- Can be operated while plugged in and recharging
- Vials can be crimped or decapped while they are in most standard removable sample trays
- Adjustable crimp settings for compatibility with most vial/septum/seal combinations including aluminum, steel and bi-metallic seals
- Crimp force sensing automatically determines when a proper seal has been formed and opens the jaws to release the vial

Cat. No.	Description	Unit
Crimpers		
70-100050	8mm Electronic Crimper 6.4 Volt	Each
70-100051	11mm Electronic Crimper 6.4 Volt	Each
70-100052	13mm Electronic Crimper 6.4 Volt	Each
70-100053	20mm Electronic Crimper 6.4 Volt	Each

Decappers		
70-100054	11mm Electronic Decapper 6.4 Volt	Each
70-100055	13mm Electronic Decapper 6.4 Volt	Each
70-100056	20mm Electronic Decapper 6.4 Volt	Each

Accessories		
70-100057	Lithium Ion Battery 6.4 Volt	Each



Greyhound Chromatography & Allied Chemicals
6 Kelvin Park, Birkenhead, Merseyside, CH41 1LT, UK
Tel: (+44) 0 151 649 4000 • Fax: (+44) 0 151 649 4001
Email: info@greyhoundchrom.com • Web: www.greyhoundchrom.com