FIRENZE - ITALIA

PHARMACY DIVISION

PRODUCTS CATALOGUE

MC-1/2 – MC-3/4

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MC-2 manual filling machines to process 100 capsules from size 000 to 5. The machines allow the simultaneous opening, filling and closing procedure for 100 capsules. Combined with manual orientator MC-1, which is also available to process capsule sizes 000 to 5, and which is capable of orienting in two phases 50 capsules at a time on the MC-2 filling machine, it can considerably reduce working times. While the capsule opening system has been tried and tested for many years, the building materials have been recently updated:

The transformation kit for changing capsule sizes is built using a special patented resin (ERTACETAL), and stainless steel for the parts in contact with the product. The frame is made of a single casting of anodized hard aluminum, so as to facilitate cleaning.

It is possible to change sizes using the transformation kit.

Moreover, upon request, accessories to facilitate the preparation are available: fixed and adjustable powder hoppers, pressure pushers of various sizes, vibrating table to facilitate the smooth descent of the powders inside the capsules (a potentiometer helps the operator to find the right intensity for the vibration).

MC3-MC-4 are, respectively, the 60-capsules-at-a-time orientator (60 + 60), and the 120-capsule filling machine, with the same characteristics described for the previous models.

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MC-2
 Production 2500 ÷ 2850 Caps./H
 Caps. Sizes 000/5
 Dimensions 18x22x10 LxWxH cm
 Mass 2 Kg
 Accessories Pressors, Hoppers, Vibrating plate

MC-3
 Production 2850 ÷ 3200 Caps./H
 Caps. Sizes 000/4
 Dimensions 19x18x7 LxWxH cm
 Mass 1 Kg
 Accessories Pressors, Hoppers, Vibrating plate

MC-4
 Production 2850 ÷ 3200 Caps./H
 Caps. Sizes 000/4
 Dimensions 18x25x10 LxWxH cm
 Mass 2 Kg
 Accessories Pressors, Hoppers, Vibrating plate

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Studied for the galenic laboratory of the Pharmacy, the MC 100 system is made up of an orientator and of a capsule filling machine mounted on the same conveniently sized base that can fit on any work table.

The capsule fitting machine is equipped with a vibrating plate and clamps to facilitate processing the powders and the full range of accessories allows to process of capsule sizes, both traditional and DB Caps.

The orientator does not require any maintenance operations and cleaning the parts of the filling machine that come in contact with the powders is extremely easy.

The MC-50 is also available upon request, with the same structure and the same characteristics as the MC-100, but intended for the small productivity of 50 capsules.

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TECHNICAL DATA

MC-50

Batch Q.ty	50 Caps.
Production	1500 Caps./H
Sizes	000-00-0-1-2-3-4-5 DB Caps AAA-AA-A-B-C-D-E
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Power Consumption	70 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	67x40x50 LxWxH cm
Mass	30 Kg

MC-100

Batch Q.ty	100 Caps.
Production	3200 Caps./H
Caps. Sizes	000-00-0-1-2-3-4-5 DB Caps AAA-AA-A-B-C-D-E
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Power Consumption	70 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	67x40x50 LxWxH cm
Mass	40 Kg

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Set of machines designed to answer the needs of the most demanding laboratories that diligently process two capsule sizes.

The system is made up of two capsule filling machines equipped with vibrating plate and pressor and with the accessories for two capsule sizes and an orientator suitable for both filling machines are installed. Said solution allows to have the filling machines always available in the different sizes, thus ensuring no time is wasted during the adjustments every time the accessories are changed. Moreover, with this set it is possible to devote a machine exclusively for one type of preparation, for example phytotherapeutic, leaving the second machine for the other normal master preparations.

The machines are positioned on single bases so as to make them easily movable on the work table or, on occasion, easily storable when not in use.

Just like the MC 100 from which they derive, the ordinary maintenance and cleaning operations are extremely easy and require very little time.

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A semi-automatic 150-capsule filling machine suitable for pharmaceutical laboratories for master preparation or for small production.

It is equipped with a vibrating plate and clamps for compressing powders.

The machine is suitable for all the sizes of hard gelatine capsules from 000 to 5 and, by using special accessories, also for DB Caps formats.

The procedure for changing format is extremely simple and quick. Some accessories with which it is possible, for example, to reduce the number of holes to produce 60 or 90 capsules or special hoppers to insert granules or pellets into capsules contribute to make the machine even more versatile.

The machine can be cleaned easily and the accessories coming into contact with the powders can also be placed in the autoclave.

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TECHNICAL DATA

Combination with	MS-4/7
Production	4000 Caps./H
Sizes	000-00-0-1-2-3-4-5 DB Caps AAA-AA-A-B-C-D-E
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Power Consumption	80 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	49x35x55 LxWxH cm
Mass	32 Kg

Combination with	MS-1
Production	6000 Caps./H
Sizes	000-00-0-1-2-3-4-5 DB Caps AAA-AA-A-B-C-D-E
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Power Consumption	80 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	49x35x55 LxWxH cm
Mass	32 Kg

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Semi-automatic capsule filling machine for preparing 300 capsules simultaneously.

Normally combined with the MS 4/7 or MS 1 orientator, the machine is suitable for all sizes of hard gelatine (from 000 to 5) and, by using special accessories, also for DB Caps formats.

The machine is equipped with a vibrating plate with adjustable intensity and a unit for compressing the powders that facilitates their smooth descent inside the capsules.

The procedure for changing format is extremely simple and quick. Some accessories with which it is possible, for example, to reduce the number of holes or special hoppers to insert granules or pellets into capsules make the machine even more versatile.

The machine can be cleaned easily and quickly and the accessories coming in contact with the powders can be removed extremely easily, thus facilitating the operators' daily maintenance.

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TECHNICAL DATA

Combination with	MS-4/7
Production	8000-10000 Cpas./H
Sizes	000-00-0-1-2-3-4-5 DB Caps AAA-AA-A-B-C-D-E
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Power Consumption	80 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	49x35x62 LxWxH cm
Mass	40 Kg

Combination with	MS-1
Hourly Production	11000-13000
Sizes	000-00-0-1-2-3-4-5 DB Caps AAA-AA-A-B-C-D-E
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Power Consumption	80 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	49x35x62 LxWxH cm
Mass	40 Kg

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The MS 4/7 capsule orienter is used to position and to evenly orientate the hard gelatine capsules on the relevant plates before they are open and filled. MS 4/7 is normally combined with capsule filling machines MS 3N, 6N and 6AUT.

A special trolley with alternate shifting makes it possible to fill both the 150-hole plates and the 300-hole ones taking up very little space on the work station.

The machine reaches the Customer equipped with accessories to process one capsule size, but by means of very simple operations it is possible to change the grids to process the different sizes from 000 to 5 and greater, in addition to the special DB Caps.

It is very easy to clean and requires no maintenance operations.

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Pneumatically-driven semi-automatic capsule filling machine to make 300 capsules.

The opening and closing operations of the capsules and the compression of powders are carried out by pressing buttons located at the front of the machine and directly connected to the pneumatic unit. This ensures that the operations run perfectly and, most importantly, that the compression of powders inside the capsules is always consistent.

Normally combined with the MS 4/7 or MS 1 orientator, the machine is available for all hard gelatine capsule sizes from 000 to 5 and, with special accessories, also for the DB Caps sizes. The procedure for changing size is extremely practical and quick and requires, after the fitting has been changed, a simple adjustment.

The cleaning operations are also very simple: the accessories coming in contact with the powders can be dislodged extremely easily, facilitating the operators' daily maintenance duties.

As optional it is possible supply silenced compressor.

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TECHNICAL DATA

Combination with	MS-4/7
Production	10000 Caps./H
Sizes	000-00-0-1-2-3-4-5
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Required Compressed Air	0,7 Mpa
Power Consumption	85 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	50x40x70 LxWxH cm
Mass	55 Kg

Combination with	MS-1
Production	16000 Caps./H
Sizes	000-00-0-1-2-3-4-5
Accessories	Capsule reducers, Hoppers to fill tablets pellets, etc
Required Compressed Air	0,7 Mpa
Power Consumption	85 Watt
Power Supply	110/230 Volt Single Phase
Dimensions	50x40x70 LxWxH cm
Mass	55 Kg

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MS-1 is an automatic orientator capable of quickly orienting 150 or 300 capsule at a time in less than a minute.

This machine is ideal used in combination with one or more semi-automatic capsule filling machines such as models MS-3N, MS-6N and MS-6AUT and is used mainly in R&D labs and by pharmaceutical, phitotherapeutic and herbal companies with medium production.

The parts coming in contact with the product can be easily inspected and dislodged, and are made of FDA approved materials. The shift of production is extremely easy and quick.

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Multigel Srl