

**CAUTION:** When filling or manually clearing the installed syringe in the PB600, do not engage the dispenser button, as the index rod will jam the syringe plunger.

After successfully filling the syringe, firmly press the dispenser button until it bottoms out, before releasing it. Repeat this procedure until the syringe is empty.

When using the optional disposable tips with LT or TLL syringes, install the optional fill stop (31) on the index rod to prevent over-filling the disposable tips, allowing sample into the syringe.

**CAUTION:** In actual use, the proper way to hold the PB600 is with the thumb on the dispenser button and the index and middle fingers on the opposite side of the body assembly. Do not operate the PB600 while holding it in the palm of the hand with the fingers wrapped around the syringe barrel. Excessive pressure on the glass barrel will fracture the syringe at the glass flange, resulting in permanent damage.

## CLEANING AND MAINTAINING THE PB600 REPEATING DISPENSER

Under normal circumstances, the PB600 Repeating Dispenser requires little or no maintenance. The dispenser should be operated in a "dry" condition; that is, without any grease, light oils, or graphite lubricants applied to the index rod or dispenser button areas. Occasionally clean the index rod with a stiff brush and acetone to keep the teeth of the index rod free of dust and dirt particles. Do not disassemble the PB600 assembly for cleaning.

If the force required to press the dispenser button needs to be modified for a specific application, the friction plug set screw (14) may be adjusted. Do not back out the set screw too far and remove all the force on the friction plug or the unit may not index correctly. Repeat step 6 of the Syringe Installation procedure after making an adjustment to verify that the PB600 is functioning properly.

## RETURN OF GOODS

Hamilton Company's return and repair policy is written to protect its employees from potentially hazardous materials (e.g., serum, radioactive materials, carcinogenic chemicals, etc.) or any substance that may cause them partial or permanent disability during the inspection or repair process. In returning a product, the customer acknowledges that the product is free from any hazardous materials. Furthermore, the customer assumes responsibility should the returned product be determined to be hazardous.

### Trademarks:

GASTIGHT is a registered trademark of Hamilton Company.  
MICROLITER is a registered trademark of Hamilton Company.

**HAMILTON**  THE MEASURE OF EXCELLENCE™

Hamilton Company manufactures products for precision fluid measuring: syringes, valves diluter/dispensers, and robotic sample processors, as well as polymeric HPLC columns and resins. For additional information any of these product lines, please contact Hamilton Company or your local Hamilton representative.

**Hamilton Company**  
4970 Energy Way Reno, Nevada 89502  
USAToll-Free 800-648-5950 Fax  
+1-775-856-7259 Telephone  
+1-775-858-3000 e-mail:  
sales@hamiltoncompany.com

**Hamilton Bonaduz AG**  
P.O. Box 26CH-  
7402 Bonaduz/Switzerland Fax  
+41-81-660-60-70 Telephone  
+41-81-660-60-60 e-mail:  
hamopd@bonaduz.hamilton.ch

<http://www.hamiltoncompany.com>

**HAMILTON**   
THE MEASURE OF EXCELLENCE™

## PB600-1 REPEATING DISPENSER

*Congratulations! You have purchased the finest quality precision fluid measuring device available today. We at Hamilton Company combine top quality materials with skilled work-manship, ensuring the highest possible performance level of every precision fluid device we manufacture. With proper care and handling, the PB600-1 Repeating Dispenser will provide unsurpassed performance in precision liquid handling year after year.*

The Hamilton PB600 Repeating Dispenser is among the most versatile, hand-held, semi-automated dispensers available. The PB600-1 dispenser has virtually unlimited applications when there is a need to deliver liquid volumes of 0.5 µL to 50 µL repeatedly up to 50 times before refilling. Some typical uses of the PB600 include preparing standard solutions, filling microtiter trays, feeding cell cultures, spotting TLC plates, dispensing VDRL antigen, human leukocyte locus-A (HLA) tissue typing and dosing small animals. Industrial applications include lubricating small bearings or parts, applying liquid solder, and dispensing mercury.

## WARRANTY STATEMENT

Hamilton Company unconditionally guarantees its products to be free of defects in materials and workmanship. Any product which fails due to such defects will be repaired or replaced at our discretion without cost, provided the device is returned with an explanation. It is the responsibility of the purchaser to determine the suitability of application and material compatibility of the products based on the published specifications of the products.

Syringes and needles manufactured by Hamilton Company are intended for scientific research and laboratory use only and are not intended for human in vivo use.

## PARTS AND ACCESSORIES

Numbers in parentheses refer to part numbers in Figure 1.

### REPLACEMENT PARTS

PB6001 .....	83700
(3) Screw 6-32 x 3/8 .....	16204
(4) Plunger Arm for 700/1700 Series Syringes .....	15144
(5) Plunger Arm for 1001/1001.25/1002 Series Syringes .....	15142
(7) PB600-1 Flange Screw .....	14704
(8) O-ring 2-12 .....	16102
(9) Index Rod .....	15104
(10) Index Rod E-ring 19661 .....	
(11) Friction Plug .....	15111

(12) Friction Plug Spring .....	53299-01
(13) PB600 Dispenser Button Spring .....	19603
(14) Friction Plug Set Screw .....	5518-06
(31) Index Rod Fill Stop (optional) .....	15125

### ACCESSORIES FOR TLL, LT, AND RN SYRINGES

100µL Disposable Tips, 1000/pk .....	75700
1 mL Disposable Tips, 500/pk .....	75701
100µL Disposable Tip Adapter .....	31330
1 mL Disposable Tip Adapter .....	31331

*Please reference Hamilton Company's Precision Fluid Measuring Catalog for additional items.*

## SYRINGE INSTALLATION AND ASSEMBLY

The PB600-1 is shipped fully assembled (without the syringe). However, the PB600-1 package contains an additional plunger arm for smaller syringe plungers.

The following assembly procedures include numbers enclosed in parentheses that correspond to specific parts of the PB600-1. The exploded view of the PB600-1 (Figure 1) shows the locations of those parts.

1. Remove the flange screw (7), turning counterclockwise, and the O-ring (8) from the PB600 assembly (23).
2. With the index rod (9) fully extended, place the glass syringe barrel into the PB600 body. With the O-ring in the groove on the flange screw, thread the flange screw into the PB600 body. Finger tighten the screw.

**CAUTION: Over-tightening the flange screw may fracture the glass flange on the syringe barrel.**

3. Select plunger arm that will fit the selected syringe (See CAUTION below) and affix it to the index rod with the screw provided (3). Plunger arm (4) is preinstalled at the factory for 700/1700 Series syringes and (5) is for 1001, 1001.25, and 1002 Series syringes.

**CAUTION: DO NOT remove the index rod from the PB600 assembly. If the plunger arm is removed, the index rod can be removed from the PB600 assembly. When the index rod comes out of the assembly, all other parts (11, 12, 13, and 22) will fall loose. Extensive reassembly is required at this point.**

4. With the index rod (9) fully extended (without the plunger assembly installed), press and release the dispenser button (22). Repeat until the index rod no longer moves (fully engaged).
5. With the index rod fully engaged, install the syringe plunger through the hole in the plunger arm and into the syringe barrel. Adjust the plunger tip to the zero scale mark silk-screened on the syringe. Tighten screws (3).

**CAUTION:** Prior to installing any GASTIGHT® syringe plunger assembly into a glass barrel, wet the PTFE plunger tip to provide lubrication between the tip and the inner bore of the syringe barrel. Failure to wet the plunger tip may cause premature tip wear or tip leakage.

6. Grasp the plunger arm, and fully retract the plunger assembly (as if you were filling the syringe). Then press the dispenser button repeatedly, counting the number of clicks, until the index rod no longer advances. If you pressed the dispenser button less than 50 times, loosen the plunger arm screw and reposition the plunger (away from zero). Retighten the plunger arm screw. Repeat the test until you achieve 50 dispenses. No other adjustments are necessary. Your PB600 is ready for use.

**NOTE: Every time you press the dispenser button, the PB600 Repeating Dispenser mechanically advances the syringe plunger 0.047" (based on a 6 cm scale length at full volume). This action advances a fluid column within the syringe 1/50 of the syringe's volume. Therefore, when setting the plunger tip to be precisely at the zero mark. What is important is that the plunger advances 50 times.**

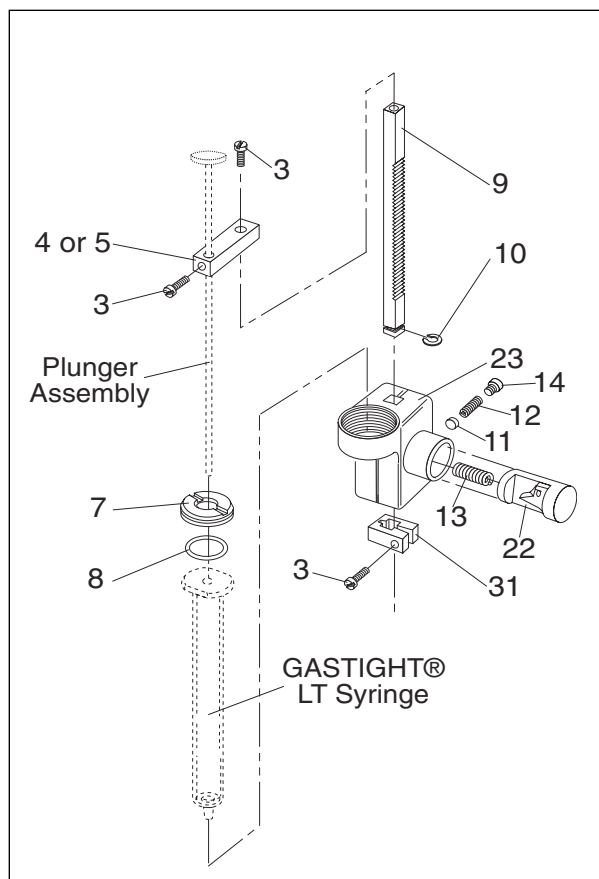


Figure 1

## SYRINGE AND NEEDLE SELECTION

Syringes and needles are not included with the PB600 Repeating Dispenser and must be ordered separately. To achieve the maximum performance from your PB600 Repeating Dispenser, consider the application, the syringe volume, and the fluid viscosity.

A syringe with a point #3 (blunt) needle is recommended for dispensing into an open vessel, into microtiter trays, or onto flat surfaces. When injecting through rubber septa, it is best to use a point #2 needle. Most fluids can be effectively handled using either a 22s-gauge or 22-gauge needle. Syringes with needles identified as (22s/2"/3) and (22/2"/3) are suitable. Hamilton part numbers for luer tip (LT), removable needle (RN), special needle (SN), standard cemented needle (N), luer tip cemented (LTN), and PTFE luer lock (TLL) syringes are identified in Table 1. Please note that LT and TLL type syringes do not come with needles. Refer to Hamilton Company's Product Catalog for more detailed information.

In addition, there are at least two special applications for the PB600-1 repeating dispenser:

**HLA Applications** - Use syringes specified in the parts list with needles designated in parentheses: (22/2"/3). For a "shooting" syringe needle information will appear as follows: (22s/2"/3). The "dropping" syringe needle appears as (22/2"/3).

**VDRL Applications** - Special luer tip syringes are designed to dispense 1/100 ml, 1/75 ml, or 1/60 ml when installed in the PB600. The syringes dispense VDRL antigen emulsion and buffered saline solutions. A special 18-gauge, point #3 (blunt) bent needle (P/N 90432) and other gauged hypodermic type needles are based on Antigen manufacturers' VDRL data sheets.

## SYRINGE AND NEEDLE SELECTION

Syringes and needles are not included with the PB600 Repeating Dispenser and must be ordered separately. To achieve the maximum performance from your PB600 Repeating Dispenser, consider the application, the syringe volume, and the fluid viscosity.

A syringe with a point #3 (blunt) needle is recommended for dispensing into an open vessel, into microtiter trays, or onto flat surfaces. When injecting through rubber septa, it is best to use a point #2 needle. Most fluids can be effectively handled using either a 22s-gauge or 22-gauge needle. Syringes with needles identified as (22s/2"/3) and (22/2"/3) are suitable. Hamilton part numbers for luer tip (LT), removable needle (RN), special needle (SN), standard cemented needle (N), luer tip cemented (LTN), and PTFE luer lock (TLL) syringes are identified in Table 1. Please note that LT and TLL type syringes do not come with needles. Refer to Hamilton Company's Product Catalog for more detailed information.

In addition, there are at least two special applications for the PB600-1 repeating dispenser:

**HLA Applications** - Use syringes specified in the parts list with needles designated in parentheses: (22/2"/3). For a "shooting" syringe needle information will appear as follows: (22s/2"/3). The "dropping" syringe needle appears as (22/2"/3).

**VDRL Applications** - Special luer tip syringes are designed to dispense 1/100 ml, 1/75 ml, or 1/60 ml when installed in the PB600. The syringes dispense VDRL antigen emulsion and buffered saline solutions. A special 18-gauge, point #3 (blunt) bent needle (P/N 90432) and other gauged hypodermic type needles are based on Antigen manufacturers' VDRL data sheets.

Table 1. Recommended Hamilton syringes for use with the PB600-1

Vol. to Dispense μL	Syringe Vol. μL	700 Series MICROLITER					1700 & 1000 Series GASTIGHT					
		N Pt. 2 22s	SN Pt. 3 22	SN Pt. 3	RN Pt. 2	LT	N Pt. 2	LTN Pt. 2	RN Pt. 2	RN Pt. 3	LT	TLL
0.5	25	80400	80465	80419	80430	80401	80200		80230	80265	80201	
1.0	50	80500	80565	80521	80530	80501	80900		80930	80965	80901	80920
2.0	100	80600	80665	80621	80630	80601	81000		81030	81065	81001	81020
5.0	250	80700	80765		80730	80701	81100		81130	81165	81101	81120
10.0	500	80800	80865		80830	80801		81217	81230	81265	81201	81220
20.0	1,000							81317	81330		81301	81320
25.0	1,250							82017			82001	
50.0	2,500							81417	81430		81401	81420
VDRL-mL												
1/100	500											81201
1/75	666.5											87801
1/60	833.5											87800

## FILLING AND DISPENSING

To avoid the formation of bubbles, slowly fill the syringe. If bubbles occur, follow one of these steps:

1. Invert the PB600 assembly and tap the glass syringe barrel to release the bubbles into the needle end of the syringe. Then press the dispenser button.
2. Manually clear the syringe by freely pushing the plunger of the syringe until the liquid is expelled.